

## Transporting, Installing, Passing On and Disposing of the Centrifuge 5430 R

These instructions for use complement the operating manual for the Centrifuge 5430 R and do not replace it. You should therefore also read the operating manual before using the centrifuge for the first time. The current version of the operating manual can be found on the Internet at [www.eppendorf.com](http://www.eppendorf.com).

### Transport



The centrifuge 5430 R may only be transported by two persons. For transportation over longer distances, a transport aid, e.g. a sack barrow, must be used.

### Installation



The weight of the centrifuge is 56 kg.  
The installation of the device must always be completed by two people.

#### 1. Open the carton.

- Cut the adhesive tape.
- Open all 4 flaps of the carton.
- Remove all accessories, rotors and packing material.



#### 2. Lift the device out of the carton.

- Lift and remove the transport securing material at the front and rear.
- Two persons should lift the centrifuge out of the carton with a firm grip on the textile straps.



#### 3. Place device on a stable, horizontal and resonance-free lab bench.



## Transporting, Installing, Passing On and Disposing of the Centrifuge 5430 R

### 4. Pull off the textile straps (do not cut) and remove the plastic covering.

- After installation, wait four hours before switching the centrifuge on in order to avoid damage to the compressor.
- Observe a minimum distance of 30 cm to adjoining devices at the sides and 15 cm from the rear side to the wall.
- Install the device in a well-ventilated location which is protected from direct sunlight to prevent it from overheating.



### 5. Check that the mains voltage and frequency match the requirements on the device type plate (see right side of the device) and then connect the device to the power supply.



### 6. Switch on the device at the mains power switch. The device is now ready to use.

### 7. Retain the packaging and all parts of the transport securing material for shipping the device at a later date.

#### Packing

Pack the centrifuge in the reverse order.

Transport of the device is only permitted in the complete and undamaged original packaging.

#### Passing On the Device

When passing the device on to third parties, be sure to include the operating manual and these instructions for use.

#### Disposal



When disposing of the device, observe the applicable legal regulations.

The disposal is regulated within the European Community by national regulations based on EU Directive 2002/96/EC on waste electrical and electronic equipment (WEEE).

For more information, please contact your supplier.

**eppendorf**  
*In touch with life*

Your local distributor: [www.eppendorf.com/worldwide](http://www.eppendorf.com/worldwide)

Eppendorf AG · 22331 Hamburg · Germany · Phone: +49 40 538 01-0 · Fax: +49 40 538 01-556 · E-mail: [eppendorf@eppendorf.com](mailto:eppendorf@eppendorf.com)

Eppendorf North America, Inc. · One Cantiague Road · P.O. Box 1019 · Westbury, N.Y. 11590-0207 · USA

Tel: +1 516 334 7500 · Toll free phone: +1 800 645 3050 · Fax: +1 516 334 7506 · E-mail: [info@eppendorf.com](mailto:info@eppendorf.com)

#### Application Support

Europe, International: Phone: +49 1803 666 789 · E-mail: [support@eppendorf.com](mailto:support@eppendorf.com)

North America: Phone: +1 800 645 3050 ext. 2258 · E-mail: [support\\_na@eppendorf.com](mailto:support_na@eppendorf.com)

Asia Pacific: Phone: +60 3 8023 6869 · E-mail: [support\\_asiapacific@eppendorf.com](mailto:support_asiapacific@eppendorf.com)